

Стерильные разъемы ASEPTIQUIK X

Технические характеристики

По вопросам продаж и поддержки обращайтесь:

Алматы (727)345-47-04
Ангарск (3955)60-70-56
Архангельск (8182)63-90-72
Астрахань (8512)99-46-04
Барнаул (3852)73-04-60
Белгород (4722)40-23-64
Благовещенск (4162)22-76-07
Брянск (4832)59-03-52
Владивосток (423)249-28-31
Владикавказ (8672)28-90-48
Владимир (4922)49-43-18
Волгоград (844)278-03-48
Вологда (8172)26-41-59
Воронеж (473)204-51-73
Екатеринбург (343)384-55-89

Иваново (4932)77-34-06
Ижевск (3412)26-03-58
Иркутск (395)279-98-46
Казань (843)206-01-48
Калининград (4012)72-03-81
Калуга (4842)92-23-67
Кемерово (3842)65-04-62
Киров (8332)68-02-04
Коломна (4966)23-41-49
Кострома (4942)77-07-48
Краснодар (861)203-40-90
Красноярск (391)204-63-61
Курск (4712)77-13-04
Курган (3522)50-90-47
Липецк (4742)52-20-81

Магнитогорск (3519)55-03-13
Москва (495)268-04-70
Мурманск (8152)59-64-93
Набережные Челны (8552)20-53-41
Нижний Новгород (831)429-08-12
Новокузнецк (3843)20-46-81
Ноябрьск (3496)41-32-12
Новосибирск (383)227-86-73
Омск (3812)21-46-40
Орел (4862)44-53-42
Оренбург (3532)37-68-04
Пенза (8412)22-31-16
Петрозаводск (8142)55-98-37
Псков (8112)59-10-37
Пермь (342)205-81-47

Ростов-на-Дону (863)308-18-15
Рязань (4912)46-61-64
Самара (846)206-03-16
Санкт-Петербург (812)309-46-40
Саратов (845)249-38-78
Севастополь (8692)22-31-93
Саранск (8342)22-96-24
Симферополь (3652)67-13-56
Смоленск (4812)29-41-54
Сочи (862)225-72-31
Ставрополь (8652)20-65-13
Сургут (3462)77-98-35
Сыктывкар (8212)25-95-17
Тамбов (4752)50-40-97
Тверь (4822)63-31-35

Тольятти (8482)63-91-07
Томск (3822)98-41-53
Тула (4872)33-79-87
Тюмень (3452)66-21-18
Ульяновск (8422)24-23-59
Улан-Удэ (3012)59-97-51
Уфа (347)229-48-12
Хабаровск (4212)92-98-04
Чебоксары (8352)28-53-07
Челябинск (351)202-03-61
Череповец (8202)49-02-64
Чита (3022)38-34-83
Якутск (4112)23-90-97
Ярославль (4852)69-52-93

Россия +7(495)268-04-70

Казахстан +(727)345-47-04

Беларусь +(375)257-127-884

Узбекистан +998(71)205-18-59

Киргизия +996(312)96-26-47

эл.почта: ccb@nt-rt.ru || сайт: <https://cpc.nt-rt.ru>

ASEPTIQUIK® X SERIES CONNECTORS

AseptiQuik® X Large Format 1" Connectors provide quick and easy sterile connections for high flow applications, even in non-sterile environments. AseptiQuik® X connectors' "TWIST-PULL-TWIST" design enables users to quickly transfer large volumes of media easily with less risk of operator error. The connectors' robust design provides reliable performance without the need for clamps, fixtures or tube welders. Biopharmaceutical manufacturers can now make 1" sterile connections with the quality and market availability they expect from the leader in single-use connection.



SPECIFICATIONS

OPERATING PRESSURE

Up to 60 psi, 4.1 bar

OPERATING TEMPERATURE

39°F to 104°F (4°C to 40°C)

STERILIZATION

Standard (blue) version:

Gamma: Up to 50 kGy irradiation

High Temperature (white) version:

Gamma: Up to 50 kGy irradiation

Autoclave: One cycle up to 266°F (130°C) for 30 minutes

TERMINATIONS

3/4" and 1" ID hose barb (19.0mm and 25.4mm), 1-1/2" sanitary

MATERIALS

Main Components:

Polycarbonate (white)

Lock Ring:

Polycarbonate, PVDF (blue)

Pull Tabs:

Polycarbonate (blue, standard version)

Polycarbonate (white, HT version)

Caps:

Polypropylene (clear)

Seals:

Silicone (clear), platinum-cured

Membrane:

Polyethylene (standard version)

Hydrophobic polyethersulfone (HT version),

PTFE strip sticker

WARNING: Pressure, temperature, chemicals, and operating environment can affect the performance of couplings. It is the customer's responsibility to test the suitability of CPC's products in their own application conditions.

FEATURES

TWIST-PULL-TWIST design

Membrane pull tabs

Robust construction

Integrated lock ring

CPC Click

BENEFITS

Intuitive three-step connection process reduces risk of operator error

Ensure simultaneous and secure removal of both membranes


Repeatable and reliable performance with no additional hardware required

Secures final connection preventing disassembly

Audible confirmation of completed assembly steps

TYPICAL FLOW RATE:

Cv Value Range: 15-55

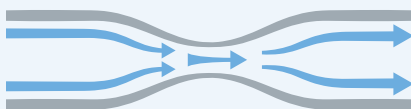

Cv values represent the approximate expected flow rate in gallons per minute of water at room temperature for a 1 PSI pressure drop. The flow is generally constrained by the smallest diameter, which in some cases will be the termination diameter and not the Nominal Flow Path.


DID YOU KNOW

Did you know that the AseptiQuik L Series Connectors also have 3/4", 1", and 1 1/2" sanitary connections?

Scan code to visit webpage


ASEPTIQUIK® X SERIES DIMENSIONS

COUPLING BODY Polycarbonate - with Blue Pull Tabs

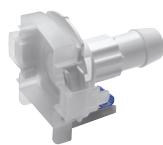

TERMINATION	METRIC EQ.	PART NO.	A	B	C
3/4" HOSE BARB	19.1 mm	AQX17012	3.54"	3.93"	1.50"
1" HOSE BARB	25.4 mm	AQX17016	3.54"	4.04"	1.53"
1-1/2" SANITARY	38.1 mm	AQX33024	3.54"	4.18"	1.73"

COUPLING BODY Polycarbonate - HT with White Pull Tabs

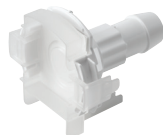

TERMINATION	METRIC EQ.	PART NO.	A	B	C
3/4" HOSE BARB	19.1 mm	AQX17012HT	3.54"	3.93"	1.50"
1" HOSE BARB	25.4 mm	AQX17016HT	3.54"	4.04"	1.53"
1-1/2" SANITARY	38.1 mm	AQX33024HT	3.54"	4.18"	1.73"

COUPLING INSERT Polycarbonate - with Blue Pull Tabs




TERMINATION	METRIC EQ.	PART NO.	A	B	C
3/4" HOSE BARB	19.1 mm	AQX22012	3.76"	4.34"	1.50"
1" HOSE BARB	25.4 mm	AQX22016	3.76"	4.45"	1.53"
1-1/2" SANITARY	38.1 mm	AQX44024	3.76"	4.58"	1.74"

COUPLING INSERT Polycarbonate - HT with White Pull Tabs




TERMINATION	METRIC EQ.	PART NO.	A	B	C
3/4" HOSE BARB	19.1 mm	AQX22012HT	3.76"	4.34"	1.50"
1" HOSE BARB	25.4 mm	AQX22016HT	3.76"	4.45"	1.53"
1-1/2" SANITARY	38.1 mm	AQX44024HT	3.76"	4.58"	1.74"

PRODUCT DIMENSIONS

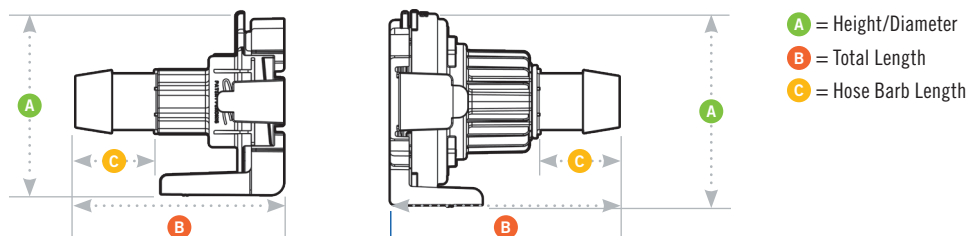

AQX ASSEMBLY PROCEDURE

STEP 1

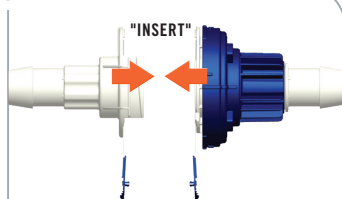


Align male and female connectors using the START alignment feature of the blue lock ring with the rib indicator of the white body. Insert the two halves together.

STEP 2



Twist the blue lock ring clockwise (approximately 75°) until audible "CPC Click" is heard. Alignment of the small blue lock ring rib indicator with the white body's rib indicator confirms final rotation is complete.

STEP 3

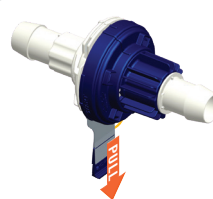

Snap the membrane pull tabs together and pull from connector.

STEP 4



Twist the blue lock ring clockwise until the final audible "CPC Click" is heard (90° from the initial starting point). Alignment of the long blue lock ring rib indicator with the white body's rib indicator confirms initial rotation is complete.

Алматы (727)345-47-04
Ангарск (3955)60-70-56
Архангельск (8182)63-90-72
Астрахань (8512)99-46-04
Барнаул (3852)73-04-60
Белгород (4722)40-23-64
Благовещенск (4162)22-76-07
Брянск (4832)59-03-52
Владивосток (423)249-28-31
Владикавказ (8672)28-90-48
Владимир (4922)49-43-18
Волгоград (844)278-03-48
Вологда (8172)26-41-59
Воронеж (473)204-51-73
Екатеринбург (343)384-55-89

Иваново (4932)77-34-06
Ижевск (3412)26-03-58
Иркутск (395)279-98-46
Казань (843)206-01-48
Калининград (4012)72-03-81
Калуга (4842)92-23-67
Кемерово (3842)65-04-62
Киров (8332)68-02-04
Коломна (4966)23-41-49
Кострома (4942)77-07-48
Краснодар (861)203-40-90
Красноярск (391)204-63-61
Курск (4712)77-13-04
Курган (3522)50-90-47
Липецк (4742)52-20-81

Магнитогорск (3519)55-03-13
Москва (495)268-04-70
Мурманск (8152)59-64-93
Набережные Челны (8552)20-53-41
Нижний Новгород (831)429-08-12
Новокузнецк (3843)20-46-81
Новокузнецк (3496)41-32-12
Новосибирск (383)227-86-73
Омск (3812)21-46-40
Орел (4862)44-53-42
Оренбург (3532)37-68-04
Пенза (8412)22-31-16
Петрозаводск (8142)55-98-37
Псков (8112)59-10-37
Пермь (342)205-81-47

Ростов-на-Дону (863)308-18-15
Рязань (4912)46-61-64
Самара (846)206-03-16
Санкт-Петербург (812)309-46-40
Саратов (845)249-38-78
Севастополь (8692)22-31-93
Саранск (8342)22-96-24
Симферополь (3652)67-13-56
Смоленск (4812)29-41-54
Сочи (862)225-72-31
Ставрополь (8652)20-65-13
Сургут (3462)77-98-35
Сыктывкар (8212)25-95-17
Тамбов (4752)50-40-97
Тверь (4822)63-31-35

Тольятти (8482)63-91-07
Томск (3822)98-41-53
Тула (4872)33-79-87
Тюмень (3452)66-21-18
Ульяновск (8422)24-23-59
Улан-Удэ (3012)59-97-51
Уфа (347)229-48-12
Хабаровск (4212)92-98-04
Чебоксары (8352)28-53-07
Челябинск (351)202-03-61
Череповец (8202)49-02-64
Чита (3022)38-34-83
Якутск (4112)23-90-97
Ярославль (4852)69-52-93

Россия +7(495)268-04-70

Казахстан +(727)345-47-04

Беларусь +(375)257-127-884

Узбекистан +998(71)205-18-59

Киргизия +996(312)96-26-47

По вопросам продаж и поддержки обращайтесь:

эл.почта: ccb@nt-rt.ru || сайт: <https://cpc.nt-rt.ru>